



3 Laboratory benches and sinks

In our new **SCALA** laboratory furniture system, laboratory benches are of major significance.

The consequent separation of services supply and furniture creates flexibility in the laboratory.

All variants of our benches can be selected with various worktop materials for a large number of application possibilities everywhere in the laboratory.

High stability, straightforward design and perfect appearance characterise our laboratory benches.

Access to water must meet various requirements in the laboratory.

Large sink modules, integrated sinks, drip cups and sink modules in service modules or fume cupboards are integrated in the laboratory as required by the specific situation.

Wherever mobility is required, our mobile units are used: under the service wing, for the suspended service boom, the service columns and the service ceiling – for fast moving in the laboratory.



Laboratory benches	108	Special tables	127
Combinations of materials/bench frames.....	108	Add-on table for low level fume cupboards.....	127
Worktop material.....	110	Balance table.....	128
Bench with H-frame.....	115	Rack.....	129
Bench with C-frame.....	116	Heavy duty rack.....	130
Bench with cantilever frame.....	117	Swing.....	131
Bench with self-supporting underbench units.....	118	Height-adjustable table.....	132
Mobile tables	119	Round table.....	133
Laboratory sinks	120	Sliding element Sekretär.....	134
Laboratory sink.....	120	Sliding element Assistent.....	135
Laboratory sink module.....	121	Sliding element Protector.....	136
Drip cup on service spine.....	123		
Drip cup in worktop.....	124		
Mobile sink.....	125		
AquaEL.....	126		



3 Laboratory benches and sinks

Our benches offer a large number of possible applications.

Our new bench frames are made of precision rectangular tubes with reinforced cross-section. The bench frames can carry a load of 200 kg without any problem. Optimally protected against external effects through the entirely homogenous powder coating, our bench frames have a flawless appearance.

The same applies to the surfaces of our work-tops. You can choose from our wide range of materials according to your requirements.

Bench frames for different needs

With their constructional designs, C-frame, H-frame and cantilever bench frames form the basis for our work benches depending on the requirement and application.

Different standard widths available

In order to be able to divide the workplaces in your laboratory to suit your needs, we offer a large number of frame widths.

Improved level compensation

Our new flush-mounted height-adjustable feet for C and H-frames offer up to 23 mm regulating distance, as an option up to 50 mm. Easy access and adjustment, for steady positioning.

Easy cleaning

The new height adjustment holds the C-frame approx. 30 mm above the floor. This makes cleaning the floor extremely easy.



H-frame

provides a high level of stability for add-on tables, mobile tables and analysis tables for working sitting or standing.

Underbench cabinets can be mobile or suspended and moved independent of modular size. Sitting niches are therefore possible anywhere.

C-frames

are extremely steady and can be loaded with 200 kg. They provide users with a large amount of knee and legroom with mobile and suspended underbench units.

Cantilever frame

provides the greatest legroom and lightest visual impact. It is fitted to service spines or directly to walls via its cantilever bracket design.

Suspended underbench units that can be moved

Our new profile enables underbench units suspended in cantilever and C-frames to be moved across frames.

Movable knee-hole cover panels

For benches without underbench units we use movable and height-adjustable knee-hole cover panels. In this way, installations routed below the rear side of the table can be hidden.

Other useful helpers

Add-on tables, Swings and round tables are autonomous objects and can be combined to form new modules as required. Our height-adjustable bench can be adjusted from 700 to 950 mm.

Our multi-talent: the rack

The rack is perfect for fitting items of equipment, AquaEl and others. The robust shelves are height-adjustable and the castors enable the fast changing of location.



3 Laboratory benches and sinks

There are no limits to using sinks and drip cups in the laboratory. With a new appearance that matches our range of laboratory furniture **SCALA** and is made of tried-and-tested materials, our sink elements can be perfectly integrated where they are needed. Materials such as stoneware, polypropylene, stainless steel and epoxy are extremely durable.

Stoneware sink modules

Our sink modules can be integrated as end sink units or along the service spine. The module made of high-strength baked and glazed stoneware in 1200 mm width is made of one piece without joints. Our sink modules are mounted on plinth units that can be fitted with drawers and hinged or tilting doors as desired.

Sink modules and drip cups

Sink modules made of stoneware or polypropylene are integrated into the service spine above the bench. Drip cups are fitted directly in the worktop. They are made of stoneware, polypropylene, epoxy resin or stainless steel.



Laboratory sink

Sinks are permanently installed components of laboratory furniture and placed against the service spine or a wall. Sinks can be combined with various types of worktop materials in many versions.






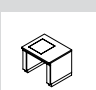



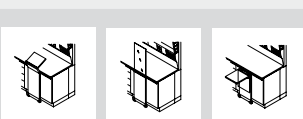
Mobile sink and AquaEl

The mobile sink with castors supplements the variable laboratory below the service wing and service ceiling. The mobile sink is connected directly to the service wing or service ceiling system via flexible pipes. AquaEl is a ready to plug in compact system for the easy supply and disposal of water in service modules. A lifting unit disposes of the waste water through the respective system.

Laboratory benches

Combinations of materials/bench frames

Combinations of worktop and bench frame materials

Material, worktop	Coated particle-board	Coated particle-board (postforming)	Solid grade laminate	Solid grade laminate Trespa Toplab+	Polypropylene
 H-frame	X	X	X	X	X
 C-frame	X	X	X	X	X
 Cantilever bench frame	X	X	X	X	X
 Mobile table frame	X	X	X	X	X
 H-frame for low level fume cupboards	-	-	-	-	X ⁴⁾
 Balance table	X	X	-	-	-
 Swing	X ¹⁾	-	-	-	-
 Round table	X ¹⁾	-	-	-	-
 Rack	X ²⁾	-	-	-	-
 Sliding elements	X ³⁾	-	-	-	-

¹⁾ Walnut veneer or light grey

²⁾ Shelves white, top of Sekretär walnut veneer

³⁾ Only walnut veneer


⁴⁾ Material with surrounding increased edge


Material, worktop	Epoxy	Stainless steel	Stoneware	Composite worktop	Glass
 H-frame	x	x	x	x	x
 C-frame	x	x	x	x	x
 Cantilever bench frame	x	x	x	x	x
 Mobile table frame	x	x	x	x	x
 H-frame for low level fume cupboards	x ⁴⁾	x ⁴⁾	x ⁴⁾	-	-


⁴⁾ Material with surrounding increased edge


Laboratory benches

Worktop material

Melamine resin facing/postforming	
Critical substances	Acids in concentrations > 10 %
Damaging substances	Concentrated hydrochloric acids Nitric acid Heated sulphuric acid
Advantage	Flat
Limitations	Joints sensitive to moisture Medium chemical resistance
Use	Mobile table, add-on table, window benches Instrument benches and laboratory benches in the dry area Cannot be used in the moist or wet area
Weight [kg/m ²]	19.6
Overall thickness [mm]	30
	Light grey NCS S 3005 R80B


Polypropylene	
Critical substances	Hydrocarbons Citric acid Oxalic acid Carbon tetrachloride Diesel oil
Damaging substances	Ozone Concentrated nitric acid Chloroform Petrol Benzol
Advantage	No joints Flat Light High chemical resistance to acids and many solvents Easy to dispose of Less breakage of glass
Limitations	Soft surface sensitive to scratches Sensitive to heat
Use	Areas with high resistance to chemicals Working with hydrofluoric acid Radio-isotope area Areas in which the lack of joints is important
Weight [kg/m ²]	20.3
Overall thickness [mm]	30
Increased edge [mm]	7
	Light grey NCS S 3005 R80B


Solid grade laminate	
Critical substances	Acids in concentrations > 10 %
Damaging substances	Concentrated hydrochloric acids Nitric acid Heated sulphuric acid
Advantage	Moisture-resistant Flat Easy to dispose of
Limitations	Reduced coating thickness
Use	Wet rooms Physical laboratories Tables with average load
Weight [kg/m ²]	26.4
Overall thickness [mm]	19
	Light grey NCS S 3005 R80B


Solid grade laminate Trespa Toplab+	
Critical substances	Acids in concentrations > 10 %
Damaging substances	Concentrated hydrochloric acids Nitric acid Heated sulphuric acid
Advantage	Antibacterial Highly-compressed surface structure High chemical resistance Moisture-resistant Flat Easy to dispose of
Limitations	Reduced coating thickness
Use	Chemical, microbiological, genetical-engineering laboratories
Weight [kg/m ²]	26.4
Overall thickness [mm]	19
	Glacier blue Similar to NCS 1010 R80B


Laboratory benches

Worktop material

Epoxy	
Critical substances	Various solvents Diluted acids
Damaging substances	Hydrofluoric acid Concentrated warm mineral acids
Advantage	No joints Flat Solid panel High mechanical load capacity Easy to dispose of
Limitations	Surface sensitive to scratches Sensitive to concentrated acids
Use	Laboratory workstation of all type
Weight [kg/m ²]	32
Overall thickness [mm]	19
Increased edge [mm]	7
	Platinum grey Similar to NCS S 4202-R

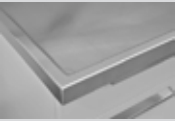
Stainless steel	
Critical substances	Cadmium Lactic acid Oxalic acid
Damaging substances	Compounds containing chlorine and bromine Formic acid Sulphuric acid
Advantage	No joints High resistance to solvents High temperature resistance
Limitations	Sensitive to halogens and their compounds
Use	For maximum loads in the area of decontamination and moisture resistance as well as solvent resistance Biology Microbiology Pharmacy Radio-isotope area Pathology
Weight [kg/m ²]	27.5
Overall thickness [mm]	30
Increased edge [mm]	7
	

Stoneware	
Critical substances	None
Damaging substances	Hydrofluoric acid
Advantage	Best chemical resistance Mechanically stable Easy to dispose of
Limitations	Evenness tolerances due to firing process Thermodynamic stress limited
Use	Areas subject to very high chemical and mechanical stress
Weight [kg/m ²]	56
Overall thickness [mm]	26
Increased edge [mm]	7
	Light grey NCS S 3005 R80B

Composite worktop	
Critical substances	None
Damaging substances	Hydrofluoric acid
Advantage	Flat Lighter than stoneware Best chemical resistance Easy to dispose of
Limitations	Thermodynamic stress limited
Use	Areas with very high chemical stress
Weight [kg/m ²]	40
Overall thickness [mm]	30
Increased edge (epoxy resin) [mm]	7
	White Similar to NCS S 0300-N

Laboratory benches

Worktop material

Glass	
Critical substances	None
Damaging substances	Hydrofluoric acid
Advantage	Flat High chemical resistance
Limitations	Sensitive to knocks at corners and edges
Use	Laboratory benches of all types subject to large amounts of chemicals
Weight [kg/m ²]	38
Overall thickness [mm]	30
	Light green NCS S 2010 G10Y


Laboratory benches

Bench with H-frame

Intended use

- Bench frame with worktop made of various materials as a work surface and storage area for laboratory work
- Supporting construction for analytical equipment and superstructures

Design

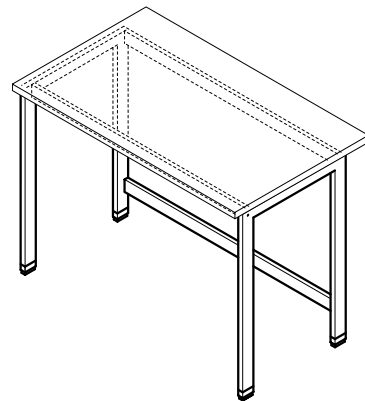

Technical data

Dimensions					
Width [mm]	600	900	1200	1500	1800
Depth [mm]	600 750 900				
Working height [mm]	750 900				

Load bearing capacity	
H-frame [kg]	200 (for fixing to the wall or for fixing to a service spine)

Design characteristics	
Construction	For suspended underbench units, cannot be moved for all kinds of frames For underbench units on castors
Height-adjustable feet	Individually adjustable

Material	
Bench frame	Steel profile 60/25/2 mm
Worktop	Depending on requirement
Height-adjustable feet	Plastic housing with steel spindle


Laboratory benches

Bench with C-frame

Intended use

- Bench frame with worktop made of various materials as a work surface and storage area for laboratory work
- Supporting construction for analytical equipment and superstructures

Design

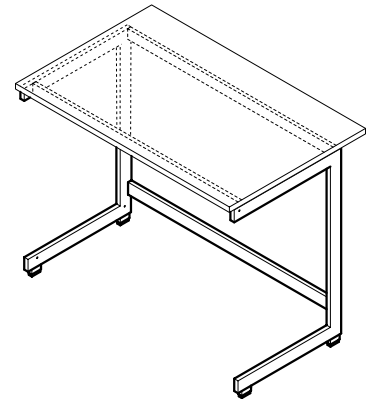

Technical data

Dimensions					
Width [mm]	600	900	1200	1500	1800
Depth [mm]				600	
				750	
				900	
Working height [mm]				750	
				900	

Load bearing capacity	
C-frame [kg]	200

Design characteristics	
Construction	For suspended underbench units, can be moved for all kinds of frames For movable underbench units
Height-adjustable feet	Individually adjustable

Material	
Bench frame	Steel profile 70/25/3 mm
Worktop	Depending on requirement
Height-adjustable feet	Plastic housing with steel spindle


Laboratory benches

Bench with cantilever frame

Intended use

- Bench frame with worktop made of various materials as a work surface and storage area for laboratory work
- Supporting construction for analytical equipment and superstructures
- For fixing to the wall or for fixing to a service spine

Design

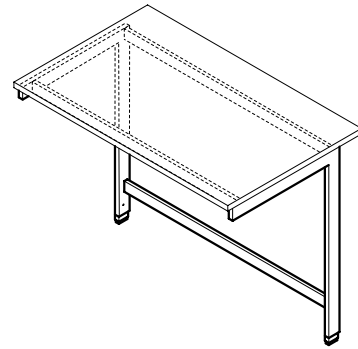

Technical data

Dimensions					
Width [mm]	600	900	1200	1500	1800
Depth [mm]			750		
			900		
Working height [mm]			750		
			900		

Load bearing capacity	
Cantilever frame [kg]	200 (for permanent mounting to a wall or a wall-mounted service spine)

Design characteristics	
Construction	For suspended and movable underbench units, can be moved for all kinds of frames
Height-adjustable feet	Individually adjustable

Material	
Bench frame	Steel profile 70/25/3 mm
Worktop	Depending on requirement
Height-adjustable feet	Plastic housing with steel spindle


Laboratory benches

Bench with self-supporting underbench units

Intended use

- Self-supporting underbench unit on plinth with worktop made of various materials as a work surface and storage area for laboratory work
- Supporting construction for analytical equipment and superstructures

Design

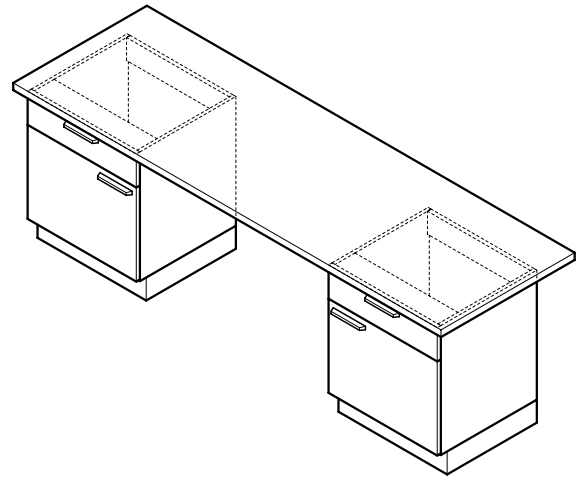

Technical data

Dimensions	
Overall width [mm]	Max. 3000
Width, underbench unit [mm]	450 600 900 1200
Total depth [mm]	750 900
Working height [mm]	750 900

Material	
Worktop	Depending on width and requirement

Load bearing capacity	
Bench with self-supporting underbench unit [kg]	200


Mobile tables

Mobile table

Intended use

- Movable bench frame with worktop made of various materials as a work surface and storage area for laboratory work
- Movable supporting construction for analytical equipment and superstructures

Design

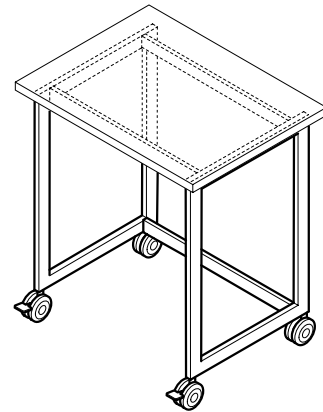

Technical data

Dimensions			
Width [mm]	900	1200	1500
Depth [mm]		600	750
		900	
Working height [mm]		750	900

Load bearing capacity	
Mobile table [kg]	150
Per heavy load castor [kg]	110

Design characteristics	
Heavy load castors	4, of which 2 can be locked (castor and steering axle can be locked)
Shelf	Optional
Underbench unit	Optional

Material	
Bench frame	Steel profile 60/25/2 mm
Worktop	Depending on requirement

Laboratory sinks


Laboratory sink

Intended use

- Water supply and disposal
- For cleaning operating equipment
- To take up large amounts of water
- Not suitable for the disposal of chemicals

Design

- 1 Underbench unit
2 Sink
3 Outlet

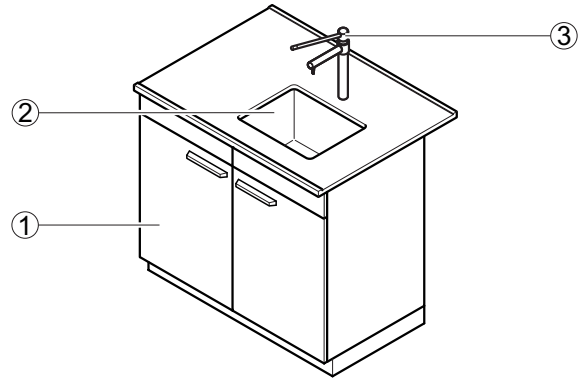

Technical data

Material Worktop	Material Sinks	Sink dimensions Width x depth x height [mm]	Type of installation
Stoneware	Stoneware	400 x 400 x 250 500 x 400 x 250	Sink installed flush with the worktop
Melamine resin facing, Solid grade laminate, Trespa Toplab+	Polypropylene	320 x 320 x 200 400 x 400 x 250 500 x 400 x 250	Sink with surrounding increased edge installed in the worktop from above
Melamine resin facing, Solid grade laminate, Trespa Toplab+	Stainless steel	340 x 370 x 150 500 x 400 x 250	Sink with surrounding increased edge installed in the worktop from above
Polypropylene	Polypropylene	385 x 385 x 250 485 x 385 x 250	Sink attached to the worktop from the bottom and welded
Stainless steel	Stainless steel	400 x 400 x 250 500 x 400 x 250	Sink welded in flush with the worktop
Composite worktop	Stoneware	380 x 380 x 250 530 x 380 x 250	Sink installed flush with the worktop
Epoxy	Epoxy	406 x 305 x 203 406 x 406 x 190 457 x 380 x 279	Sink installed flush with the worktop

Dimensions					
Width [mm]	600	900	1200	1500	1800
Depth [mm]			600 ¹⁾ 675 ¹⁾ 705 ¹⁾ 750 825 855 900		
Working height [mm]			900		

¹⁾ Positioning of the outlets on the side of the sink, if required

Sanitary technology	
Water connection	Permanent connection
Waste water connection	Permanent connection with siphon
Water fitting (tap)	Bench-mounted service outlet as an option
Eye shower	Optional


Laboratory sinks

Laboratory sink module

Intended use


- Water supply and disposal
- For cleaning operating equipment
- To take up large amounts of water
- For installation on special underbench units
- Not suitable for the disposal of chemicals

Design

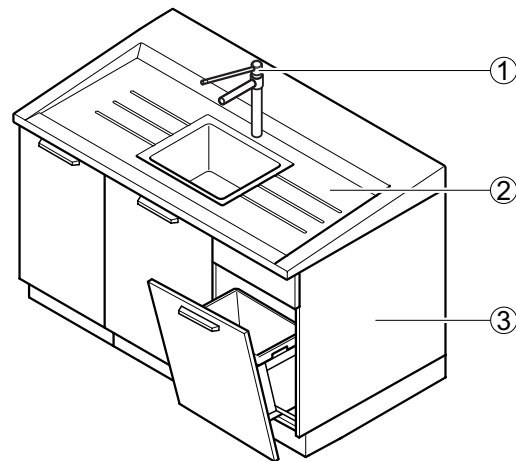


- 1 Outlet
- 2 Sink module
- 3 Underbench unit (3-piece)


Variants

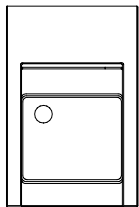


600

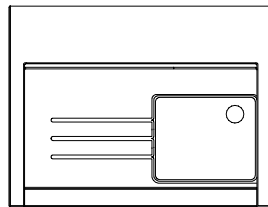


1200

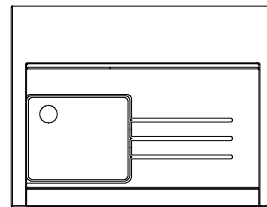

1200

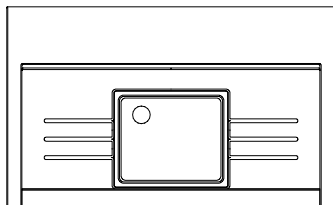

1500

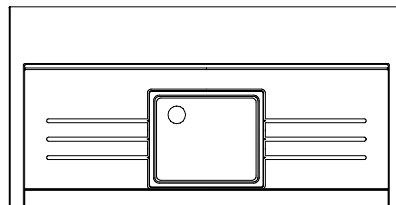

1800

Laboratory sinks

Laboratory sink module

Technical data

Dimensions				
Width, sink module [mm]	600	1200	1500	1800
Depth, sink module for wall bench with service spine [mm]	675 or 825			
Depth, sink module for double bench with service spine [mm]	705 or 855			
Depth, sink module in front of building wall [mm]	760 or 910			
Depth, sink module as an end sink [mm]	-		740	
Overall height, sink module with underbench unit [mm]	910 front to 950 rear			
Sink dimensions, width x depth x height [mm]	460 x 390 x 250			
Height, edge of sink [mm]	20 front to 50 rear			

Material

Sink module	Stoneware
-------------	-----------

Design characteristics

Construction	Self-supporting moulded draining area Surrounding increased edge
Modular design	Different underbench units possible As end sink with special underbench unit

Sanitary technology

Water connection	Permanent connection
Waste water connection	Permanent connection with siphon
Water fitting (tap)	Bench-mounted service outlets as an option
Eye shower	Optional


Laboratory sinks

Drip cup on service spine

Intended use

- Water supply and disposal
- For cleaning operating equipment
- Sink module underneath water fittings to take up small amounts of water
- Not suitable for the disposal of chemicals

Design

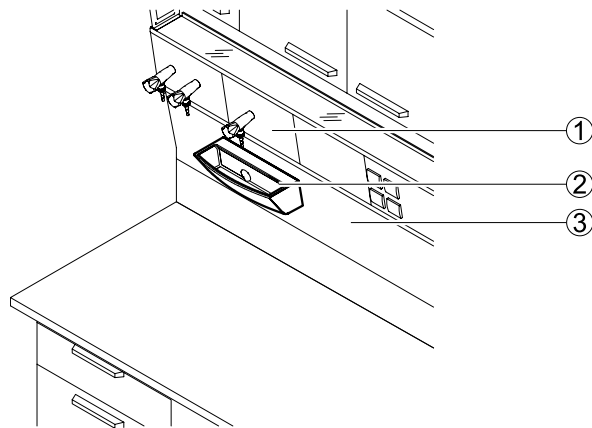

- 1 Service panel with corner valves
- 2 Sink module
- 3 Fascia panel for service spine

Technical data

Dimensions	
Width [mm]	294
Depth [mm]	132
Height [mm]	112
Internal sink dimensions width x depth x height [mm]	Approx. 270 x 85 x 80

Material	
Sink module	Stoneware Polypropylene

Design characteristics	
Construction	Attached to the fascia panel of the service spine

Sanitary technology	
Water connection	Permanent connection
Waste water connection	Permanent connection with siphon
Water fitting (tap)	Cell outlets as an option


Laboratory sinks

Drip cup in worktop

Intended use

- Water supply and disposal
- For cleaning operating equipment
- Drip cup underneath water fittings to take up small amounts of water
- Not suitable for the disposal of chemicals

Design

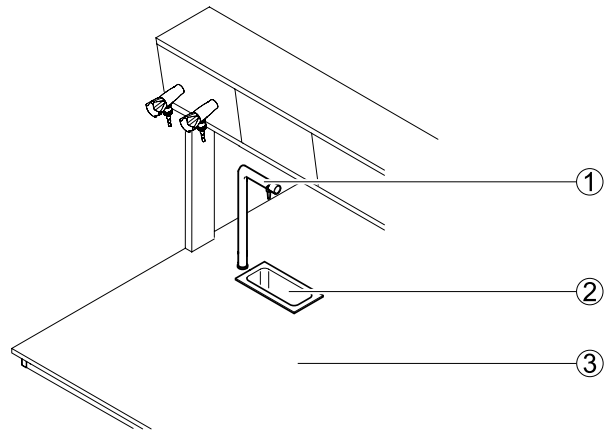

- 1 Outlet
2 Drip cup
3 Worktop

Technical data

Dimensions	
Width x depth [mm]	295 x 145
Height [mm]	Approx. 125 to 140 depending on material
Internal sink dimensions width x depth x height [mm]	Approx. 250 x 100 x 150

Material, drip cup	Material, worktop
Stoneware	Stoneware, composite worktop
Polypropylene	Polypropylene, melamine resin facing, solid grade laminate, Trespa Toplab+
Stainless steel	Stainless steel, melamine resin facing, solid grade laminate, Trespa Toplab+
Epoxy	Epoxy

Design characteristics	
Construction	Installed in the worktop from the top or bottom

Sanitary technology	
Water connection	Permanent connection
Waste water connection	Permanent connection with siphon
Water fitting (tap)	Bench-mounted service outlets as an option

Laboratory sinks


Mobile sink

Intended use

- Mobile water and gas supply and disposal
- For cleaning operating equipment at any location
- Not suitable for the disposal of chemicals

Design

- 1 Connecting pipes
- 2 Fitting with two cold water outlet points
- 3 Worktop
- 4 Sink
- 5 Underbench unit on castors

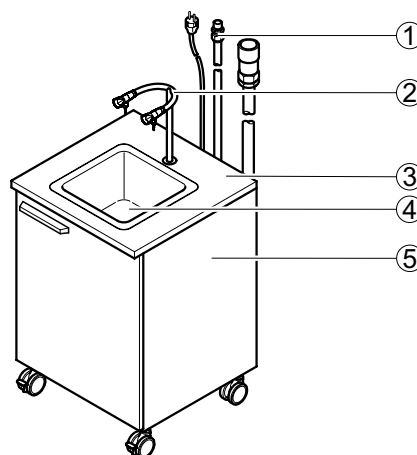

Technical data

Dimensions	
Width [mm]	605
Depth [mm]	600
Height without outlet [mm]	900
Sink dimensions width x depth x height [mm]	320 x 320 x 200
Height, [mm] castors	110
Length, supply and drain pipes [mm]	2500
Length, connecting pipes [mm]	2500

Material	
Worktop	Particle-board with melamine resin facing
Sink	Polypropylene

Load bearing capacity	
Mobile sink [kg]	150

Design characteristics	
Construction	Mounted on underbench unit on castors with hinged door Sink installed in the worktop from above Pipes and cables routed out at the rear of the underbench unit Waste water lifting unit in the underbench unit Water supply is switched off in the case of a power failure

Electrics	
Power supply [V]	230

Sanitary technology	
Water connection	Flexible with plug connector
Waste water connection	Flexible with plug connector
Gas connection	Flexible with plug connector as an option
Water fitting (tap)	Standard outlet
Gas outlet	Standard outlet combined with water fitting as an option
Mixer tap	Additional flexible water connection as an option

Laboratory sinks


AquaEL

Intended use

- Mobile water and gas supply and disposal
- For cleaning operating equipment at the workplace at any mobile or stationary laboratory workstation
- Not suitable for the disposal of chemicals

Design

- 1 Connecting pipes
- 2 Outlet with water outlet point
- 3 Housing with pump
- 4 Sink

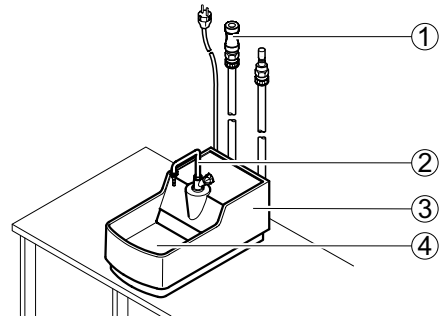

Technical data

Dimensions	
Width x depth x height (without outlet) [mm]	317 x 585 x 268
Sink, width x depth x height [mm]	260 x 275 x 105
Length, supply and drain pipes [mm]	1500
Length, connecting pipes [mm]	1500

Weight	
Weight without outlet [kg]	14

Material	
Material	GFK varnished

Design characteristics	
Construction	Compact system with flexible pipes and cables ready for connection Waste water lifting unit integrated in the housing Water supply is switched off in the case of a power failure

Electrics	
Power supply [V]	230

Sanitary technology	
Water connection	Flexible with plug connector
Waste water connection	Flexible with plug connector
Gas connection	Flexible with plug connector as an option
Water fitting (tap)	Standard outlet
Gas outlet	Standard outlet combined with water fitting as an option
Mixer tap	Additional flexible water connection as an option


Special tables

Add-on table for low level fume cupboards

Intended use

- For adding to low level fume cupboards
- Bench frame with worktop made of various materials as a work surface and storage area for laboratory work
- Supporting construction for analytical equipment and superstructures

Design

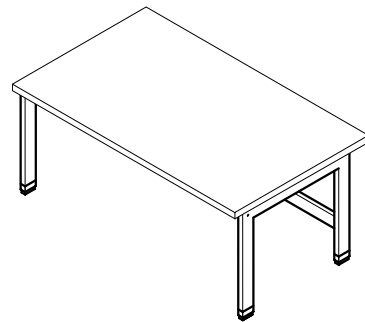

Technical data

Dimensions					
Width [mm]	900	1200	1500	1800	2100
Depth [mm]	600				
Working height [mm]	500				

Material	
Bench frame	Steel profile 60/25/2 mm
Worktop	Depending on requirement
Height-adjustable feet	Plastic housing with steel spindle

Load bearing capacity	
H-frame [kg]	200

Design characteristics	
Worktop	Surrounding increased edge
Height-adjustable feet	Individually adjustable


Special tables

Balance table

Intended use

- For setting up analytical balances and other sensitive measuring equipment
- Bench frame with worktop and specially mounted, vibration-free plate

Design

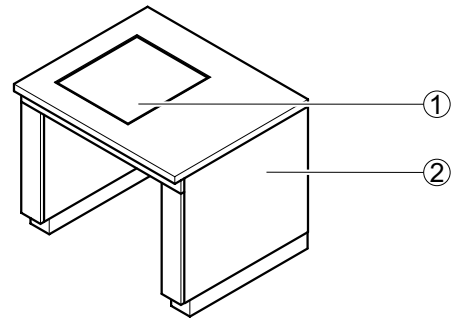

- 1 Balance plate made of fine concrete
2 Table cover

Technical data

Dimensions	
Width [mm]	900
Depth [mm]	750 900
Working height [mm]	750 900
Width x depth [mm] balance plate	400 x 450

Material	
Supporting construction	Steel profile
Worktop	Depending on requirement
Balance plate	Fine concrete


Weight	
Total weight [kg]	120
Balance plate [kg]	65

Design characteristics	
Construction	Specially mounted, heavy balance plate made of fine concrete Supporting construction with balance plate, vibration-decoupled

Intended use

- Mobile flexible storage area
- Can be used with the 600 mm deep shelf as a mobile workplace for desk work
- Not suitable for storing hazardous substances

Design

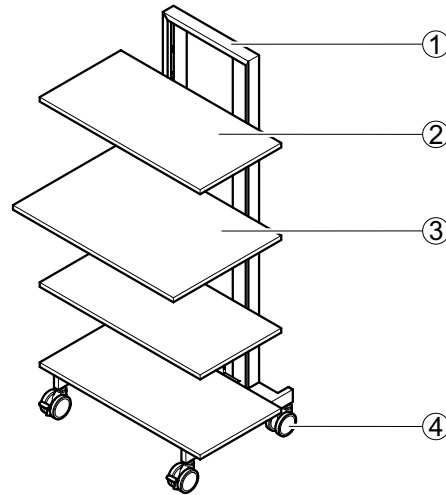

- 1 Steel support frame with grid
- 2 Shelf, depth 450 mm
- 3 Shelf, depth 600 mm
- 4 Heavy load castors with brakes

Technical data

Dimensions	
Width [mm] with shelf	900
Depth [mm] with shelf depth 450 mm	600
Height [mm]	1790
Depth, shelf [mm]	450 600

Material	
Supporting construction	Steel profile
Shelf 22 mm	Shelf of OSB board

Load bearing capacity	
Total [kg]	150
Shelf [kg]	20

Design characteristics	
Heavy load castors	4, of which 2 can be locked (castor and steering axle can be locked)
Shelf	Can be adjusted without tools with a pitch of 150 mm
Integrated distribution pillar	Optional


Special tables

Heavy duty rack

Intended use

- Mobile multi-stage storage area
- With flexible work surfaces for free horizontal configuration
- Suitable for accommodating stackable and non-stackable measuring instruments / measuring instrument towers
- Suitable for heavy apparatus

Design

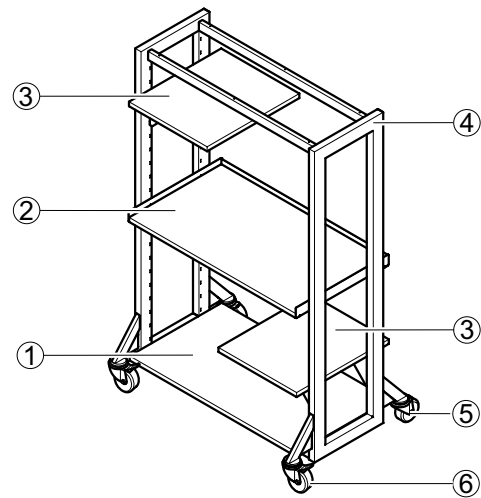

- 1 Lower shelf, fixed
- 2 Height-adjustable shelf, full width
- 3 Height-adjustable shelf, depth 590 mm
- 4 Steel support frame
- 5 Heavy load castors without brake
- 6 Heavy load castors with brake

Technical data

Dimensions		
Width [mm]	1200	1800
Depth [mm]	770	
Height [mm]	1790	
Shelf, width x depth [mm]	500 x 590 1072 x 590	500 x 590 1672 x 590

Material	
Supporting construction	Steel profile 70 x 40 mm
Shelf	Shelf of OSB board


Load bearing capacity	
Total	500 [kg]
Shelf 500 x 590 [mm]	30 [kg]
Shelf 1072 x 590 [mm]	70 [kg]
Shelf 1672 x 590 [mm]	70 [kg]
Lower shelf 1072 x 590 [mm]	150 [kg]
Lower shelf 1672 x 590 [mm]	150 [kg]

Design characteristics	
Heavy load castors	4, of which 2 can be locked (castor and steering axle can be locked)
Shelf	Can be adjusted with a grid of 75 mm

Intended use

- Bench frame with worktop made of various materials as a work surface and storage area for laboratory work
- Supporting construction for analytical equipment and superstructures
- Visually appealing connection of laboratory benches of different depths

Design

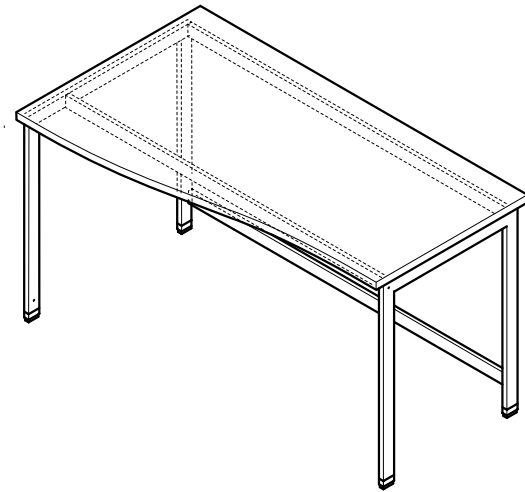

Technical data

Dimensions		
Width [mm]	1200	1500
Depth [mm]	600 – 750 750 – 900	
Working height [mm]	750 900	

Material	
Bench frame	Steel profile 60/25/2 mm
Worktop	Depending on requirement
Height-adjustable feet	Plastic housing with steel spindle

Load bearing capacity	
Total [kg]	200


Special tables

Height-adjustable table

Intended use

- Bench frame with worktop made of various materials as a height-adjustable work surface and storage area for laboratory work
- Supporting construction for analytical equipment and superstructures

Design

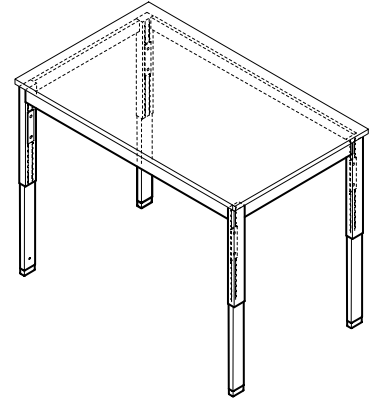

Technical data

Dimensions		
Width [mm]	1200	1500
Depth [mm]	600	750
	900	
Working height [mm]	700 – 950	

Material	
Bench frame	Steel profile 60/25/2 mm
Worktop	Depending on requirement
Height-adjustable feet	Plastic housing with steel spindle

Load bearing capacity	
Total [kg]	200

Design characteristics	
Working height	Can be adjusted with a grid of 25 mm
Bench frame	H-frame


Special tables

Round table

Intended use

- For adding to benches with H-frames and C-frames as an additional work surface

Design

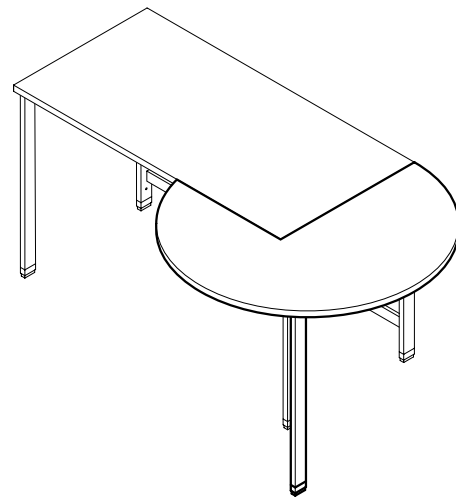

Technical data

Dimensions	
Diameter [mm]	1200
Working height [mm]	750 900

Material	
Bench frame	Steel profile 60/25/2 mm
Worktop	Depending on requirement
Height-adjustable feet	Plastic housing with steel spindle

Load bearing capacity	
Round table [kg]	50


Special tables

Sliding element Sekretär

Intended use

- Inclined desk that can be moved on a laboratory work bench

Design

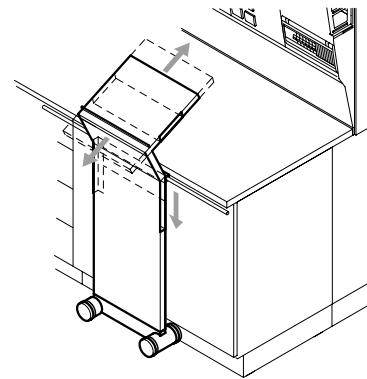

Technical data

Dimensions	
Width [mm]	416
Working height, bench [mm]	900
Max. overall height [mm]	1244


Design characteristics	
Construction	Sliding element on 4 wheels Fastened to a sliding rail on the laboratory work bench with an adjustable writing pad

Material	
Sliding element	Particle-board with melamine resin facing in walnut veneer

Intended use

- Fold-out storage area and desk that can be moved on a laboratory work bench

Design

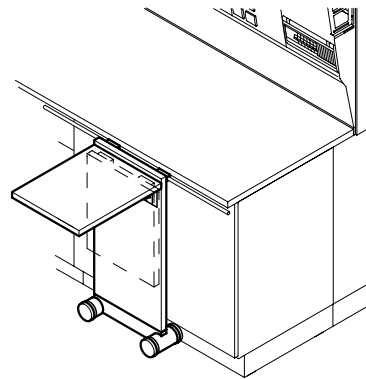

Technical data

Dimensions	
Width [mm]	406
Depth [mm]	530
Working height, bench [mm]	900

Load bearing capacity	
Storage area and desk [kg]	10

Design characteristics	
Construction	Sliding element on 4 wheels Fastened to a sliding rail on the laboratory work bench Can be folded down completely

Material	
Sliding element	Particle-board with melamine resin facing in walnut veneer


Special tables

Sliding element Protector

Intended use

- Protection from splashes and splinters that can be moved on a laboratory work bench

Design

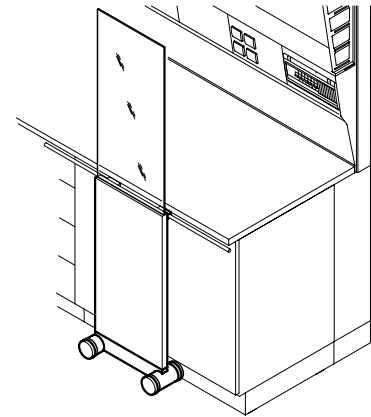

Technical data

Dimensions	
Width [mm]	406
Working height, bench [mm]	900
Overall height [mm]	1780

Design characteristics	
Construction	Sliding element on 4 wheels Fastened to a sliding rail on the laboratory work bench

Material	
Sliding element	Particle-board with melamine resin facing in walnut veneer
Splash protection	Safety glass

Laboratory benches and sinks